



QVF® is a trademark of
De Dietrich Process Systems

ROTARY EVAPORATOR

The most universal solution
for gentle and efficient evaporation






QVF® ROTADEST for ACHEMA 2018

HIGHLIGHTS

- Most universal type of evaporator
- Gentle evaporation due to film evaporation
- Evaporation down to the dry product
- Reducing foam generation due to reduced bubbling
- Avoiding hazardous bumping boiling
- Reliable industrial design

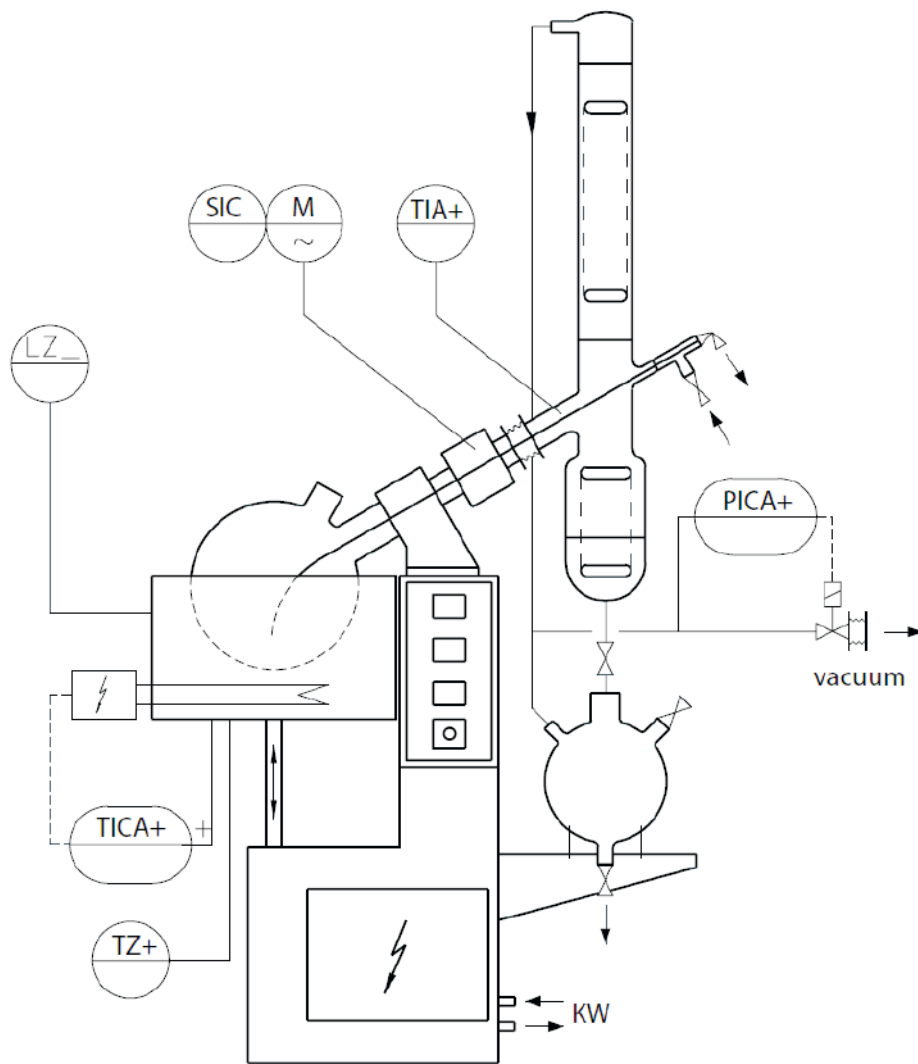

CONCEPT

Rotary evaporators are mainly used to evaporate light boiling components under normal pressure or vacuum. When used in pilot plant and small scale production they are also suitable for semi-continuous evaporation. They are characterized by providing a gentle thermal treatment for temperature-sensitive mediums and provide a higher evaporation rate as non-stirred glass distillation flasks. Compared to other evaporation processes the unique process feature of rotary evaporators is their ability to evaporate all volatile media providing at the end a dry product in the flask.

The process conditions are limited by the maximum temperature of the heating bath (max. ~ 160°C) , the temperature of the cooling liquid (e.g. ~ 15°C tap water) for the condenser and the vacuum generator (e.g.~10mbar).

The QVF® ROTADEST can be supplied for EX-rated areas up to zone 1 IIB.

FUNCTION

The liquid to be processed in the rotary evaporator is either charged batch wise by gravity, by a pump or sucked-in by vacuum into the evaporation flask. During the evaporation process further feed can be added by means of a pump, gravity or by the operating vacuum through the PTFE-hose passing the rotating shaft and ending in the spherical vessel. The rotating sphere is half way rotating through the liquid in a heating bath filled with water or another suitable liquid heating fluid. The bath is heated by means of electrical heating elements and can be lowered to immediately stop the evaporation process.

The rotation of the sphere which is flanged to the geared motor ensures an effective mixing of the product inside the vessel and wets the complete inner surface of the sphere, increasing the evaporation surface, therefore increasing the evaporation rate and hence shortening the evaporation time.

The components with the lower boiling point are evaporated in the rotating sphere first. Their vapours flow through a sleeve shaft integrated in the drive and enter the condenser assembly. The vapours are condensed and optionally cooled in a distillate cooler before being captured in a final distillate receiver.

TECHNICAL DESCRIPTION

Heating bath



The 50l and 100l rotary evaporators are supplied with a height adjustable heating bath. This heating bath is raised and lowered by a hydraulic cylinder which is operated by means of a hand pump and a pressure relief valve. As a safety feature the bath will be automatically lowered via a solenoid valve in case of a power failure or shutdown. The heating bath is heated electrically. Either water or other thermal oils can be used as a bath liquid. A transparent splash protection keeps back the bath liquid entrained by the rotation of the sphere and prevents at the same time the operator from accidentally touching the optional handhole nozzle of the rotating sphere.

Rotating spherical flask

The rotating spherical flask of 50 or 100l is made of corrosion-resistant material borosilicate glass 3.3 permitting an excellent observation of the evaporation process. The level inside the rotating flask can be monitored with a RADAR-sensor. To further intensify the mixing and for powdery substances the rotary sphere can also be supplied with inwards raised indentations. The sphere is connected to the drive by means of flap screws in order to facilitate their removal. The glass spheres are optionally equipped with a hand hole with a quick release cover to enable the addition of products directly into the sphere while being fixed to the drive.

Drive

The glass sphere of the QVF® ROTADEST is mounted to a PTFE-lined steel shaft which is connected to a gear box directly driving the rotating shaft. The gear box permits the stepless adjustment of the number of revolutions of the sphere and can be actuated by an electric motor. Optionally and alternatively an electric motor with frequency converter can be provided. The rotating PTFE-sheeted vapour shaft is sealed by a mechanical PTFE-seal at the upper end permitting the reliable self-adjusting operation for slight overpressure or low vacuum. All rotating parts including the connection flange of the rotary sphere are protected by a safety cover.

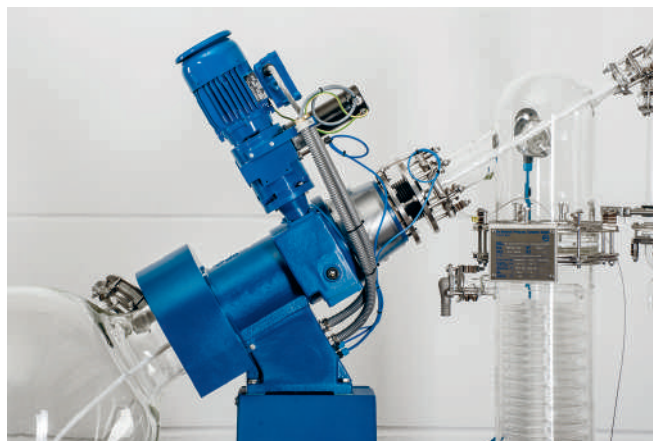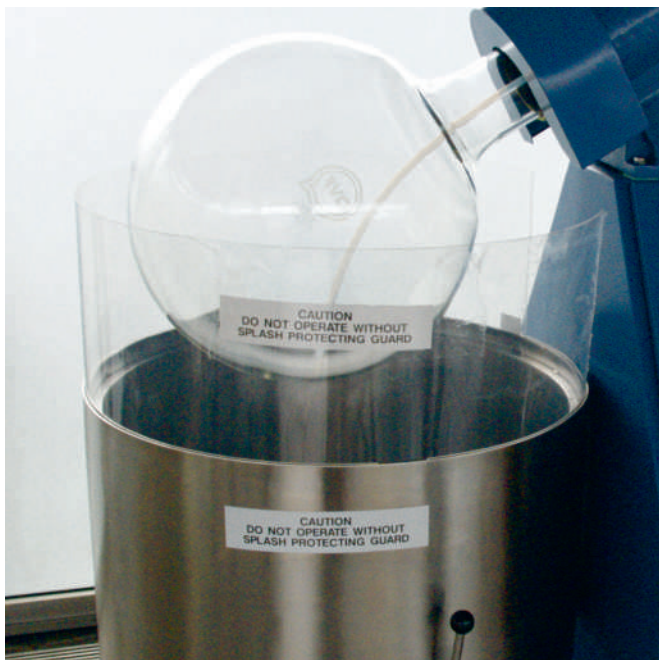


Drive with electric actuator for the gear box to adjust the speed

Condensation section

This module is the complete condenser part in which the condensate is condensed and collected. It essentially consists of a condenser with or without an after cooler, one or two recipients, the connecting pipes, including the necessary fixing as well as a vacuum connection. The condenser part is supported by a console attached to the housing. Condenser parts have already been designed and delivered in a variety of different models to cater to the special desires of the user. However, experience has shown that a wide field of application is covered by the varieties described below. They all have product feeding, concentrate and distillation extraction, connections to cooling water and vacuum as well as a ventilation and pressure measurement connection.


Condenser Version A

This condenser arrangement with integral spherical receiver has single vacuum, venting and drain connections fitted with valves. The vacuum connections is equipped with a bellows via tee-piece for connection to both vacuum pump and pressure gauge. There is also a facility for filling and suction emptying the rotating flask. This arrangement is supported by means of a bracket fixed to the cabinet. Rotary film evaporators incorporating this type of condenser arrangement are suitable for continuous operation at atmospheric pressure or batch operation under vacuum.

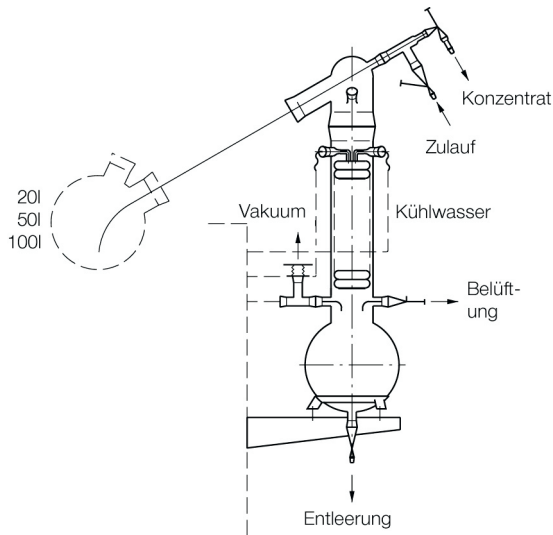


Condenser Version B

This version is basically the same as option 1 but utilises a standard condenser. This system is fitted with two spherical receivers, therefore allowing one to be exchanged, enabling continuous operation under vacuum as well as atmospheric pressure. The vacuum connection in this system is fitted with bellows to the unit.

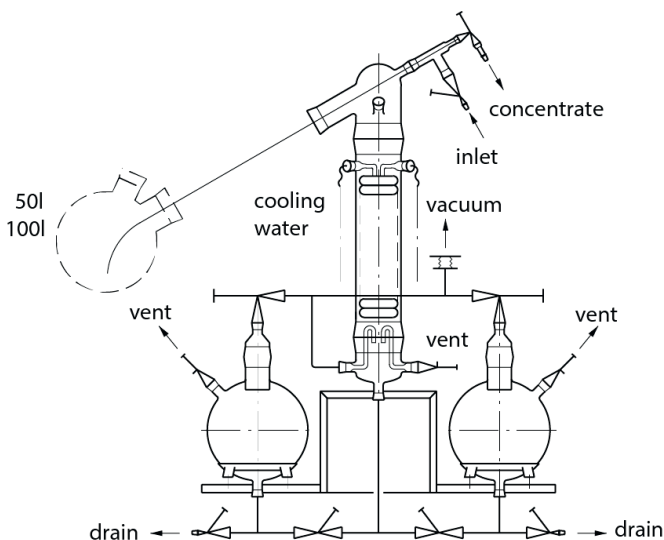

Condenser Version C

This system incorporates a uprising condenser and product cooler for particularly high cooling performance and continuous operation under vacuum. The alternately drainable receivers permit continuous operation under vacuum as well as under atmospheric pressure.

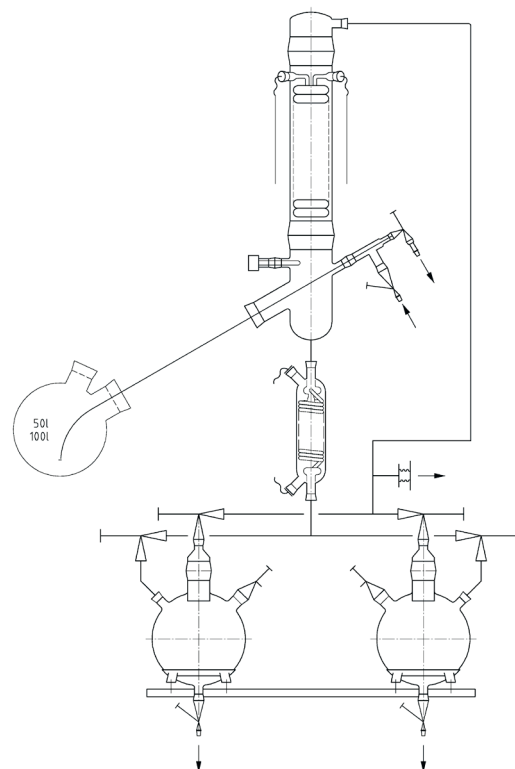

MEASURING AND CONTROL EQUIPMENT

The rotary evaporators are equipped with all necessary instrumentation to enable easy and safe operation. The operating touch panel is positioned for the convenience of the operator on one side of the upper housing part and permits full control over the process including data logging. All electronic equipment consists of standardized industrial components ensuring long term serviceability. The service access is eased from outside by a removable cover. Explosion-protected versions have an EX-proof sealed EX-housing inside the cabinet which itself is designed splash-proof (IP 54 type of protection).

Indicated and/or controlled are the following parameters:


- Head temperature in the condenser with Pt100
- Bath temperature with Pt100
- Vacuum with pressure transmitter (option)
- Liquid level inside the rotating flask via RADAR (option)
- Speed of the rotary sphere with hand wheel (remote control as option)

One power supply 230/400 V, 3 Ph, N, PE.

The QVF® ROTADEST can be installed in EX-rated areas up to zone 1 IIB.



QVF® ROTADEST with visible EX-housing for demonstration

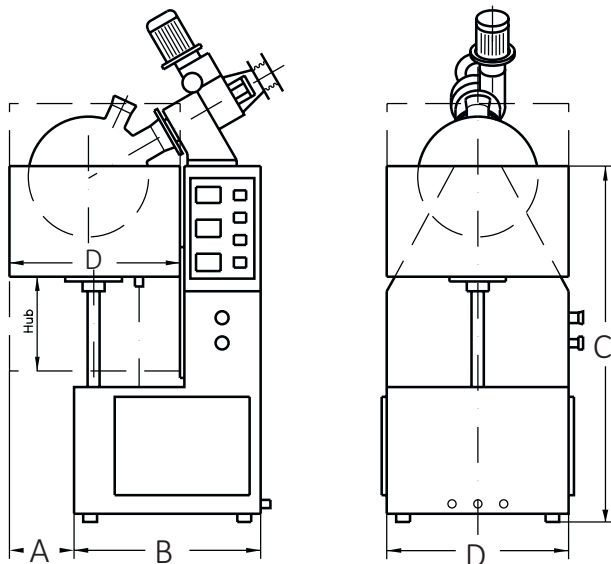


Dimensions

	Unit	RD50	RD100
A	mm	475	525
B	mm	760	810
C	mm	1450	1550
D	mm	740	840
Hub	mm	320	380


Technical Data ROTADEST

	Unit	RD50	RD100
Kolbenvolumen	l	50	100
Rotationsgeschwindigkeit	U/min	0-82	0-82
Antriebsleistung	kw	0,37	0,37
Heizleistung	kw	9	12

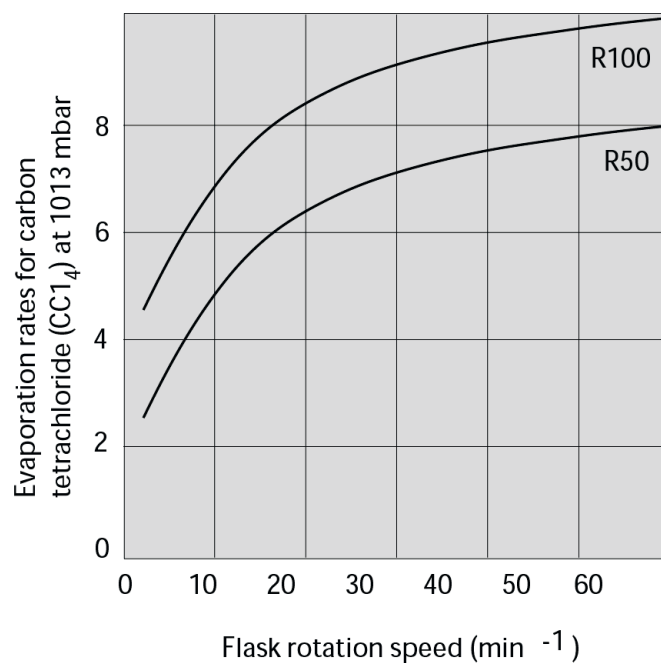

SPECIAL BENEFITS OF THE QVF® ROTADEST

- Low vacuum due to high performance mechanical PTFE-seal and certified glass flange system
- Self-adjusting mechanical PTFE-seal for slight overpressure or low vacuum
- Exclusive use of corrosion resistant materials as borosilicate glass 3.3 and PTFE for all components being in contact with the product
- FDA-Materials certificates for surfaces in contact with product
- Robust drive unit for the rotary sphere
- Condensation section suited to the process
- Proper observability of the processes
- Reliable and sustainable industrial construction
- Simple operation
- Ex-Protected version available

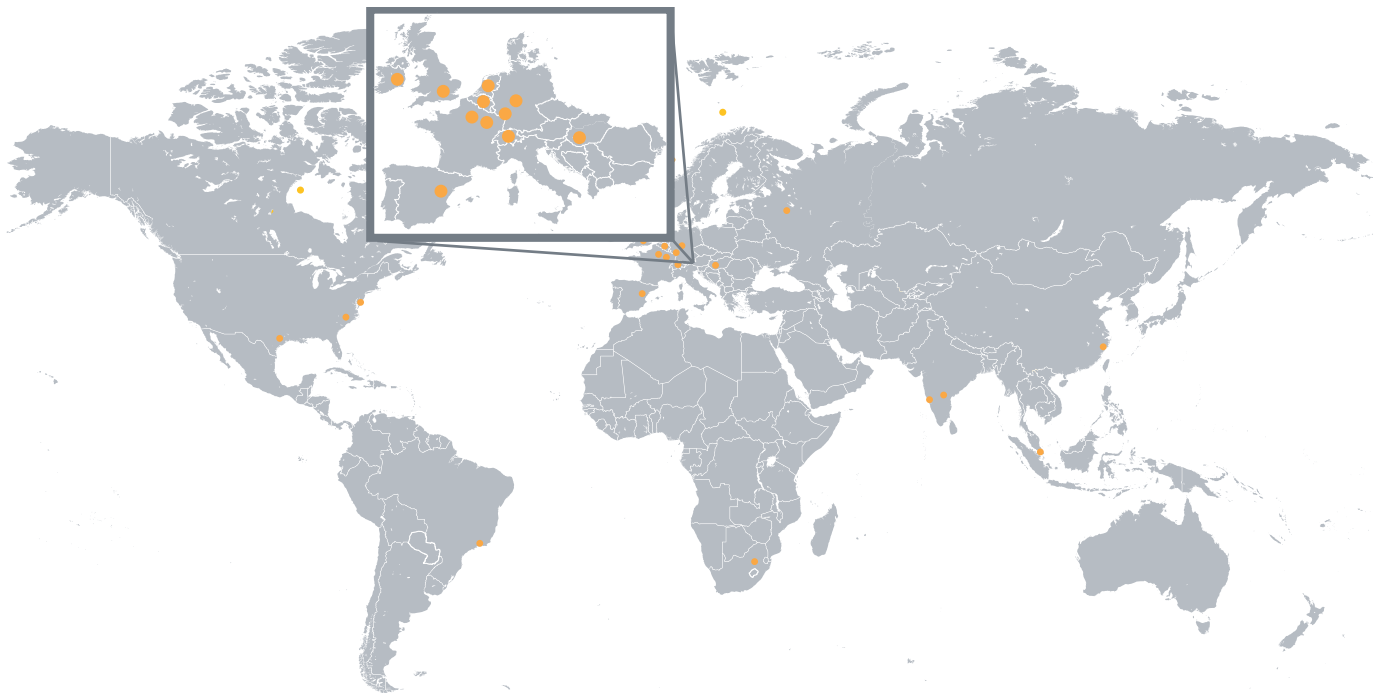

De Dietrich

PROCESS SYSTEMS



DE DIETRICH SAS

Château de Reichshoffen
F 67891 Niederbronn Cedex
Tel. +33 3 88 80 26 00
Fax +33 3 88 80 26 95
www.dedietrich.com

BENELUXSTAATEN

De Dietrich Process Systems N.V.
B - Heverlee-Leuven
Tel. +32 16 40 5000
Fax +32 16 40 5500
info@benelux.dedietrich.com

BRASILLEN

De Dietrich Do Brasil Ltda
São Paulo
Tel. +55 11 2703 7380
Fax +55 11 2702 4284
brasil@dedietrich.com.br

CHINA

De Dietrich Process Systems Co. Ltd
Wuxi
Tel. +86 510 6696 7500
Fax +86 510 6696 7599
info@dedietrichchina.com

DEUTSCHLAND

De Dietrich Process Systems GmbH
Mainz
Tel. +49 6131 9704 0
Fax +49 6131 9704 500
mail@qvf.de

FRANKREICH

De Dietrich S.A.S.
Zinswiller
Tel. +33 3 88 53 23 00
Fax +33 3 88 53 23 99
sales@dedietrich.com

De Dietrich Process Systems Semur
S.A.S.
Semur-en-Auxois
Tel. +33 3 80 97 12 23
Fax +33 3 80 97 07 58
info.semur@dedietrich.com

GROSSBRITANIEN / IRLAND

De Dietrich Process Systems Ltd
Stafford
Tel. +44 1785 609 900
Fax +44 1785 609 899
sales@ddpsltd.co.uk

INDIEN

De Dietrich Process Systems (India)
Pvt. Ltd
Mumbai
Tel. +91 22 6742 42 70
Fax +91 22 28 505 731
sales@dedietrich.co.in

Hyderabad
Tel. +91 40 2980 5442
Fax +91 40 2980 5441
sales@dedietrich.co.in

SCHWEIZ

De Dietrich Process Systems AG
Bubendorf
Tel. +41 61 925 11 11
Fax +41 61 921 99 40
info@rosenmund.com

SINGAPUR

De Dietrich Singapore (PTE) Ltd
Singapore
Tel. +65 68 61 12 32
Fax +65 68 61 61 12
info.sg@dedietrich.com

SPANIEN

De Dietrich Equipos Quimicos S.L.
Barcelona
Tel. +34 93 292 0520
Fax +34 93 21 84 709
comercial@dedietrich.es

SÜDAFRIKA

De Dietrich South Africa (PTY) Ltd
Dunswart
Tel. +27 11 918 4131
Fax +27 11 918 4133
info.za@dedietrich.com

VEREINIGTE STAATEN

De Dietrich Process Systems Inc.
Mountainside, NJ
Tel. +1 908 317 2585
Fax +1 908 889 4960
sales@ddpsinc.com

Charlotte, NC
Tel. +1 704 587 04 40
Fax +1 704 588 68 66
rosenmund@ddpsinc.com

UNGARN

De Dietrich S.A.S.
Budapest
Tel. +36 20 56 83 444
info.hu@dedietrich.com

Die internationale Unternehmensgruppe De Dietrich Process Systems ist der führende Anbieter von Systemlösungen und Reaktoren für korrosive Anwendungen sowie Anlagen für die mechanische Fest/Flüssigtrennung und Trocknung. Die Systemlösungen von De Dietrich Process Systems werden in den Industriebereichen Pharma, Chemie und Lebensmittel eingesetzt.

www.dedietrich.com